



## Shipping Configuration for the CapNMR™ probe

First, ensure your probe has been cleaned and dried out. The probe should not be shipped with solvent in the flow path. Rinse the probe with the last known solvent used, then blow out with air and then rinse with acetone and blow out with air again. Send the probe with the filters and unions attached. Use caution when restraining the capillary with the restraint wheel making sure the capillary cannot be crushed in the edge of the case.



Place the case in the shipped box (36" x 12" x 12") and insert the Styrofoam inserts along all sides of the probe case. Ensure the probe cannot shift in the box.

Place the final support sheet on top of the case and seal with common tape. Finally, make sure you have called MRM for an RGA number that you write on the shipping label. (RGA ##### – Return goods authorization followed by a 4 digit number). The shipping address will be provided when you receive your RGA number from your MRM contact person. If you have any questions, call (217) 351-4359.

